

Updating ESI SC Software (TN110610)

The following instructions may be used to update the ESI SC software to the latest version. Administrator privileges will be required on the computer.

- 1) Download software installer from: www.icpms.com/support
Login is not required to use the link at bottom of page.
- 2) Uninstall old version. (Uninstall will not remove any created FAST method files and these file will be automatically updated to the format of the new version.)
If any customization has been performed on the ESI SC software such as custom racks or special device parameters, please collect copies of these files for reference by ESI technical support.
- 3) Reboot computer.
- 4) Install from downloaded files (you may have to unzip to a folder to be able to run the files.)
 - i. If asked "Confirm Uninstall" of PL-2303 choose, no.
 - ii. If asked to "Modify? Repair? Remove?" of PL-2303 choose, "Repair".
- 5) Check in the Control Panel>>System>>Hardware>>Device Manager to verify that the computer sees any Prolific USB-Serial drivers if used. If the computer does not see them, remove them for a few seconds then plug back in. (The driver version for the Prolific Driver should be 2.0.13.130 or later).
- 6) Run ESI SC and verify that ports are configured correctly. Please refer to install instructions pertaining to the specific "ESI + host instrument" configuration.
- 7) Verify that the correct instrument/emulation is selected in ESI SC.
- 8) Place the correct racks in ESI SC layout
- 9) Initialize ESI SC and use the Manual window to verify proper operation.
- 10) Test control from host instrument software.